


DOCUMENT NUMBER: FM-000-0004-48	TITLE: Form - Product Change Notice	
REVISION: H DCO# DC0005391	DOCUMENT CLASSIFICATION: 04-Forms, Templates or Other	

PCN Number: PCN388239-0001 **Revision:** A **Creation Date:** 26 Jun 2023

This Product Change Notice supersedes the following documents:

N/A

Description and Reason for Change:

We are writing to inform you that as part of our proactive sustaining engineering program and to address component obsolescence issues, we are replacing the camera module on IntelliXcap™ Automated Screw Cap Decapper/Recapper Acoustic 96-format.

This change will be effective immediately and will affect all IntelliXcap™ Automated Screw Cap Decapper/Recapper Acoustic 96-format with Serial Numbers greater than 46-8014-1021-0012.


Impact of Change:

This change will be effective immediately and will affect all IntelliXcap™ Automated Screw Cap Decapper/Recapper Acoustic 96-format with Serial Numbers greater than 46-8014-1021-0012.

While there is no impact on image quality nor equipment performance, there is a change on the installation procedure of Decoding Software V14 used to operate the Verification Camera module (IntelliXcheck module).

Please note that this change will not impact service or support of the previous versions of the Verification camera modules.

Confidential: The information is confidential and is to be used only in connection with matters authorized by Azenta and no part of it is to be disclosed to others without prior written permission from Azenta.

DOCUMENT NUMBER: FM-000-0004-48	TITLE: Form - Product Change Notice	
REVISION: H DCO# DC0005391	DOCUMENT CLASSIFICATION: 04-Forms, Templates or Other	

Affected Part Numbers:

This PCN affects Part Numbers listed below:

Part Number	Item Description
46-8014	ASSY,INTELLIXCAP,DECAPPER,96-CHANNEL,ACOUSTIC CP 

Replacement/Alternative Part Numbers

No suitable replacement available.

Specification of alternative products

N/A.

Recommended Action:

Please follow the new installation instructions provided with IntelliXcap™ Automated Screw Cap Decapper/Recapper Acoustic 96-format when installing Decoding Software V14 on your PC.


The new installation procedure is described on document 382851 (Revision B; supplied together with the decapper).

Moreover, the patch file **uEye_cr_patch.reg** needed for camera recognition and used during the installation procedure is also supplied with the decapper and included in the provided USB stick PN 384789.


Timeframe for Implementation:

Effective immediately.

Confidential: The information is confidential and is to be used only in connection with matters authorized by Azenta and no part of it is to be disclosed to others without prior written permission from Azenta.

DOCUMENT NUMBER: FM-000-0004-48	TITLE: Form - Product Change Notice	
REVISION: H DCO# DC0005391	DOCUMENT CLASSIFICATION: 04-Forms, Templates or Other	

Appendix:

	Installation of Azenta Decoding Software and Setup of Verification Camera	382851 Revision B
---	---	----------------------

Technical Information

Electrical Category	Type 2 - Equipment is energized. Energized circuits are covered or insulated
Materials	USB stick (PN: 384789)

Safety Requirements

NOTICE

Untrained or Improperly Equipped Personnel

- Untrained or improperly equipped personnel performing this procedure may cause damage to the equipment.
- Personnel performing this procedure must read and understand this procedure and have the proper tools and supplies ready before starting.
 - Personnel performing this procedure must know the applicable safety codes, facility safety procedures, safety equipment, and emergency contact information.

Installation


Step	Action
1.	Power ON your PC and connect the provided USB stick to your PC. <i>NOTE: Make sure your computer meets the minimum system requirements of the program:</i> <i>Processor: Intel Core i5, 8GB RAM</i> <i>Free Disk Space: minimum of 1GB</i>
2.	Ensure you have full admin rights (Read, Write, and Modify) to the registry. Close all other applications during the installation process. <i>NOTE: Do not connect the IntelliXcap for Acoustic to the PC at this stage.</i>
3.	Open your file navigator and navigate to the connected USB drive.
4.	Right-click on <code>azenta_decoding_software_setup_14_xx</code> and select Run as administrator .
5.	Click Next on the Welcome screen to start the install process.

Proprietary information - This document and the information disclosed herein is confidential and proprietary to Azenta, Inc. and may not be reproduced in whole or in part or disclosed to any third party or used without the prior written consent of Azenta, Inc.

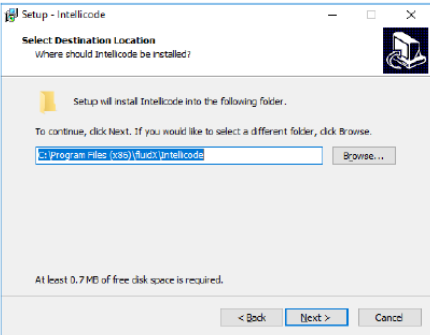
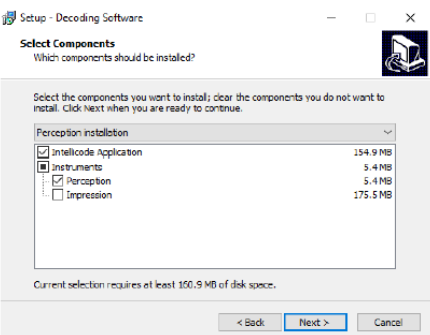
www.azenta.com
Service and Sales Support: +1.888.2.AZENTA

Page Number 1 of 5
Part Number 382851 Rev. B

Confidential: The information is confidential and is to be used only in connection with matters authorized by Azenta and no part of it is to be disclosed to others without prior written permission from Azenta.

DOCUMENT NUMBER: FM-000-0004-48	TITLE: Form - Product Change Notice	
REVISION: H DCO# DC0005391	DOCUMENT CLASSIFICATION: 04-Forms, Templates or Other	

Setup and Configuration Procedure
IntelliXcap Automated Acoustic Cap Decapper


Step	Action
6.	<p>Verify the install path and click Next. The installation proceeds.</p>  <p>The Select Components window appears.</p>
7.	<p>Ensure the IntelliCode Application and Perception check boxes are selected and then click Next.</p> 

Proprietary information - This document and the information disclosed herein is confidential and proprietary to Azenta, Inc. and may not be reproduced in whole or in part or disclosed to any third party or used without the prior written consent of Azenta, Inc.

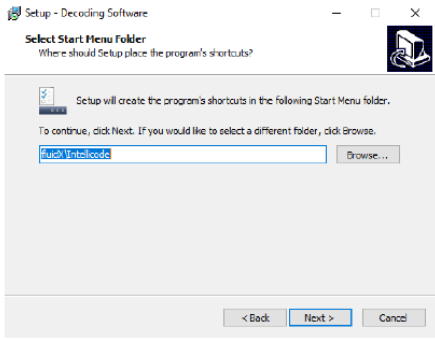
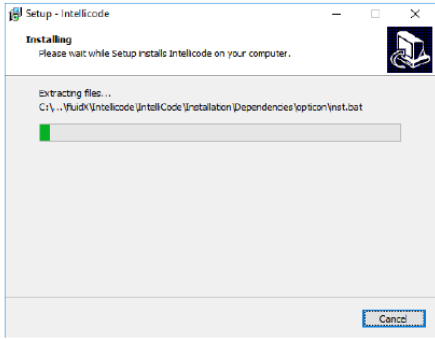
www.azenta.com
Service and Sales Support: +1.888.2.AZENTA

Page Number 2 of 5
Part Number 382851 Rev. B

Confidential: The information is confidential and is to be used only in connection with matters authorized by Azenta and no part of it is to be disclosed to others without prior written permission from Azenta.

DOCUMENT NUMBER: FM-000-0004-48	TITLE: Form - Product Change Notice	
REVISION: H DCO# DC0005391	DOCUMENT CLASSIFICATION: 04-Forms, Templates or Other	

Setup and Configuration Procedure
IntelliXcap Automated Acoustic Cap Decapper


Step	Action
8.	<p>Select a shortcut location and then press Next.</p> 
9.	<p>Verify the install details and click Install. The installation proceeds.</p>  <p>The installation proceeds. This may take a few minutes.</p>

Proprietary information - This document and the information disclosed herein is confidential and proprietary to Azenta, Inc. and may not be reproduced in whole or in part or disclosed to any third party or used without the prior written consent of Azenta, Inc.

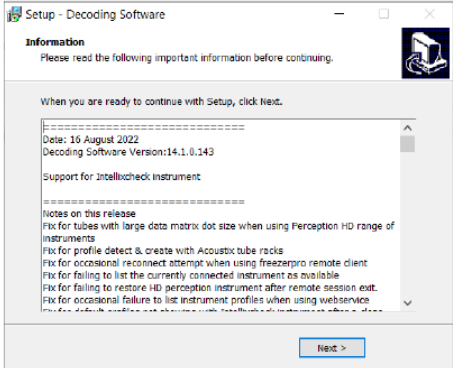
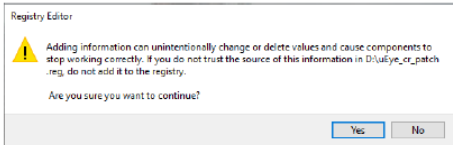
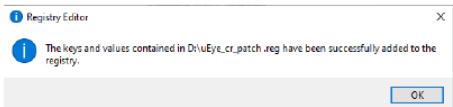
www.azenta.com
Service and Sales Support: +1.888.2.AZENTA

Page Number 3 of 5
Part Number 382851 Rev. B

Confidential: The information is confidential and is to be used only in connection with matters authorized by Azenta and no part of it is to be disclosed to others without prior written permission from Azenta.

DOCUMENT NUMBER: FM-000-0004-48	TITLE: Form - Product Change Notice	
REVISION: H DCO# DC0005391	DOCUMENT CLASSIFICATION: O4-Forms, Templates or Other	

Setup and Configuration Procedure
IntelliXcap Automated Acoustic Cap Decapper


Step	Action
10.	<p>Read the release note information and click Next.</p> 
11.	To finish the installation, select the <i>Yes, restart the computer now</i> option, and then click Finish .
12.	Confirm the Azenta Decoding Software is successfully installed on the PC.
13.	Open File Explorer and navigate to the connected USB drive.
14.	<p>Open the Decoding Software folder and right-click on the <code>uEye_cr_patch.reg</code> file, then select Merge.</p> <p>The following message is displayed.</p> 
15.	<p>Click Yes.</p> <p>The following message is displayed.</p> 

Proprietary information - This document and the information disclosed herein is confidential and proprietary to Azenta, Inc. and may not be reproduced in whole or in part or disclosed to any third party or used without the prior written consent of Azenta, Inc.

www.azenta.com
Service and Sales Support: +1.888.2.AZENTA

Page Number 4 of 5
Part Number 382851 Rev. B

Confidential: The information is confidential and is to be used only in connection with matters authorized by Azenta and no part of it is to be disclosed to others without prior written permission from Azenta.

DOCUMENT NUMBER: FM-000-0004-48	TITLE: Form - Product Change Notice	
REVISION: H DCO# DC0005391	DOCUMENT CLASSIFICATION: 04-Forms, Templates or Other	

Setup and Configuration Procedure
IntelliXcap Automated Acoustic Cap Decapper

Step	Action
16.	Click OK.
17.	Power on and plug the IntelliXcap for Acoustic device into an available port on the PC using the provided cable and wait for the device to be recognized.
18.	Open the Decoding Software. The Instrument should now be detected in the software application).

Revision History


Revision	Date
Revision A	06 JUN 2022
Revision B	21 JUN 2023

Proprietary information - This document and the information disclosed herein is confidential and proprietary to Azenta, Inc. and may not be reproduced in whole or in part or disclosed to any third party or used without the prior written consent of Azenta, Inc.

www.azenta.com
Service and Sales Support: +1.888.2.AZENTA

Page Number 5 of 5
Part Number 382851 Rev. B

Confidential: The information is confidential and is to be used only in connection with matters authorized by Azenta and no part of it is to be disclosed to others without prior written permission from Azenta.

DOCUMENT NUMBER: FM-000-0004-48	TITLE: Form - Product Change Notice	
REVISION: H DCO# DC0005391	DOCUMENT CLASSIFICATION: 04-Forms, Templates or Other	

For more information, please contact:

Sofia Peniche Maria **Senior Global Product Manager**
Tel: +44 (0) 781 596 0529 Email: sofia.peniche@azenta.com

Azenta Service and Sales Support: +1.800.379.7221.

Further contact information can be found at <https://www.azenta.com/contact-us>

1 RECORD RETENTION

For Control and Retention of Quality Records, see QMS101517

Confidential: The information is confidential and is to be used only in connection with matters authorized by Azenta and no part of it is to be disclosed to others without prior written permission from Azenta.