



# Freezer Uncrating

The reusable, quick-snapping crate reduces time, waste, and cleanup. The empty crate can be flat packed and efficiently stored or shipped back to be reused. Always inspect the freezer crate for damage and verify the bill of lading prior to accepting the shipment.

1. Locate the clip remover tool stored underneath one of the clips.
2. To remove, cover the clip with your hand and carefully rotate the tool towards the center of the panel to pry loose both the clip and the tool.
3. To remove remaining clips, insert the tool into the clip's raised pocket, cover the clip with your hand, and carefully rotate the tool towards the center of the panel.

4. Start with the top panel and proceed with the side panels.
5. Place the reinforced side panel ramp so the locating holes align with those on the skid, and then insert the mounting pins.
6. Unbolt the wheel blocks from the caster bracket and then reattach the caster bracket nuts.
7. Unlock the wheel brakes and carefully roll the freezer down the ramp.
8. Flat pack the crate and store or ship back to be reused.

